

How to Clean the Incubator

As a courtesy to the teachers and classrooms that will use the incubator after you, please clean the incubator to the best of your ability.

Responsible students can help with steps 1 and 3.

1. Discard all eggshells, unhatched eggs, water cups and paper towels.
2. With a soft sponge, use water and bleach mixture to thoroughly clean the bottom of incubator, plastic pan and hardware cloth. Rinse with hot soapy water and air dry.
3. For the lid of the incubator and the turner, use antibacterial wipes to clean off any residue and to disinfect.

